

Manual

TransforMax EPI300 Electrocompetent *E. coli* and TransforMax EPI300 Chemically Competent *E. coli*

For Research Use Only. Not for use in diagnostic procedures.

TransforMax™ EPI300™ Electrocompetent *E. coli* and TransforMax™ EPI300™ Chemically Competent *E. coli* are part of the Epicentre™ product line, known for its unique genomics kits, enzymes, and reagents which offer high quality and reliable performance.

Manual

TransforMax EPI300 Electrocompetent and Chemically Competent *E. coli*

Contents

1. Product description	3
2. Product specifications	3
3. Product designations and kit components	4
4. Storage conditions	4
5. Electroporation of TransforMax EPI300 Electrocompetent <i>E. coli</i>	4
6. Transformation of TransforMax EPI300 Chemically Competent <i>E. coli</i>	5
7. Induction of CopyControl fosmid clones	5
8. References	6
9. Technical support and product guarantee	6

Manual

TransforMax EPI300 Electrocompetent and Chemically Competent *E. coli*

1. Product description

TransforMax EPI300 *E. coli* have been specifically engineered for use with LGC, Biosearch Technologies' CopyControl™ Cloning Systems.* The cells contain a mutant *trfA* gene, whose protein product is required for initiation of replication from the *oriV* origin of replication contained on the CopyControl vectors. The *trfA* gene is under tightly regulated control of an inducible promoter. When grown on standard LB plates or in LB or SOC culture medium, expression of the *trfA* gene is repressed. Addition of the CopyControl Induction Solution (provided with the cells) induces expression of the *trfA* gene and subsequent utilisation of the *oriV* origin of replication and high copy amplification of the CopyControl fosmid clones.

Important phenotypes and applications

- Mutant *trfA* gene under tightly regulated control of an inducible promoter for complete copy number control of CopyControl clones.
- Supports blue/white screening of vectors expressing the *LacZ'* α -complementing peptide.
- Restriction minus for efficient cloning of methylated genomic (e.g. mammalian genomic) DNA.
- Endonuclease minus (*endA1*) to ensure high yields of plasmid clones.
- Recombination minus (*recA1*) to ensure the stability of large cloned inserts.

2. Product specifications

	TransforMax EPI300 Electrocompetent <i>E. coli</i>	TransforMax EPI300 Chemically Competent <i>E. coli</i>
Transformation efficiency:	Greater than 1×10^{10} cfu/ μ g DNA using 10 pg of pUC19 and an Eppendorf Multiporator (2.5 kV, fast charge rate) using a 2 mm cuvette	Greater than 5×10^8 cfu/ μ g DNA using 10 pg of pUC19 and an Eppendorf Multiporator (2.5 kV, fast charge rate) using a 2 mm cuvette.
Genotype	F ⁻ <i>mcrA</i> Δ (<i>mrr-hsdRMS-mcrBC</i>) ϕ 80d <i>lacZ</i> Δ M15 Δ <i>lacX74</i> <i>recA1</i> <i>endA1</i> <i>araD139</i> Δ (<i>ara, leu</i>)7697 <i>galU galK</i> λ ⁻ <i>rpsL nupG trfA dhfr</i>	
Quality control:	Both electrocompetent and chemically competent cells are tested for induced expression of the <i>trfA</i> gene using transformants harboring a plasmid with an <i>oriV</i> origin of replication, and to be free of contaminating DNA rendering resistance to ampicillin, tetracycline, kanamycin and chloramphenicol.	

Manual

TransforMax EPI300 Electrocompetent and Chemically Competent *E. coli*

3. Product designations and kit components

Product	Kit size	Catalogue number	Reagent description	Part number	Volume
TransforMax EPI300 Electrocompetent <i>E. coli</i>	10 x 100 µL**	EC300110	TransforMax EPI300 Electrocompetent <i>E. coli</i>	SS001009D	10 x 100 µL
			pUC19 DNA [†]	SS000200-D	10 µL
			CopyControl Induction Solution [‡]	SS000023-DCON	
	50 x 100 µL**	EC300150	TransforMax EPI300 Electrocompetent <i>E. coli</i>	SS001004-D	50 x 100 µL
			pUC19 DNA [†]	SS000200-D	10 µL
			CopyControl Induction Solution [‡]	SS000023-DCON	
TransforMax EPI300 Chemically Competent <i>E. coli</i>	10 x 50 µL**	C300C105	TransforMax EPI300 Chemically Competent <i>E. coli</i>	SS001004-D	10 x 50 µL
			pUC19 DNA [†]	SS000200-D	10 µL
			CopyControl Induction Solution [‡]	SS000023-DCON	

** Number of electroporations: EC300110: 10; EC300150: 50; C300C105: 10. EC300150 is shipped as 5 x EC300110.

[†] Each is supplied with 10 µL (100 pg/µL) of pUC19 Control DNA in TE Buffer.

[‡] CopyControl Induction Solution is also available separately as Catalogue #CCIS125 (25 mL)

4. Storage conditions

Store TransforMax EPI300 *E. coli* cells at -70 °C. The CopyControl Induction Solution and the pUC19 Control DNA can be stored at either -20 °C or -70 °C. Warm the CopyControl Induction Solution to room temperature and mix thoroughly before use.

5. Electroporation of TransforMax EPI300 Electrocompetent *E. coli*

NOTE: The electroporation procedure described here uses 50 µL of electrocompetent cells. A different volume of cells can also be used based on the experiences and needs of the user.

1. DNA should be in water or very low salt buffer (e.g. TE Buffer: 10 mM Tris-HCl [pH 7.5], 1 mM EDTA) to prevent arcing during electroporation. The pUC19 Control DNA is provided in TE at 100 pg/µL. If running a transformation control, dilute the pUC19 Control DNA 1:10 (to a final concentration of 10 pg/µL) with sterile, deionised water and use 1 µL for electroporation.
2. Prepare 1 mL of SOC or LB medium (do not include antibiotic in the medium) for each electroporation to be performed. This medium will be used for post-electroporation outgrowth of transformed cells. Maintain the medium at room temperature.
3. Pre-chill electroporation cuvettes and 1.5 mL tubes on ice.
4. Set-up the electroporation device according to the manufacturer's recommendations for bacterial (*E. coli*) electroporation.

Manual

TransforMax EPI300 Electrocompetent and Chemically Competent *E. coli*

5. Thaw TransforMax EPI300 Electrocompetent *E. coli* cells on ice on ice. Mix gently. Use the cells immediately. Unused cells can be refrozen at -70 °C.
NOTE: *Refrozen cells may have reduced transformation efficiency, but for most cases, the reduction is not significant enough to interfere with the desired results.*
6. Transfer the desired amount of DNA and 50 µL of cells to a pre-chilled microcentrifuge tube.
NOTE: *A smaller volume of cells can be used based on the needs and experiences of the user. Mix the cells and DNA by pipetting up and down 2-3 times.*
7. Transfer the cell/DNA mix to the electroporation cuvette. Be sure that there are no air bubbles in the cuvette. Wipe the cuvette of any condensation. Place into the electroporator and apply the electric pulse at the manufacturer's recommendations for bacterial (*E. coli*) electroporation.
8. Immediately after electroporation, add 950 µL of room temperature SOC medium¹ to the cuvette. Mix gently by pipetting up and down 2-3 times.
9. Transfer the cells to a 15 ml tube and incubate at 37 °C with shaking at 220-230 rpm for one hour to recover the cells and allow expression of the antibiotic resistance marker.
10. Dilute and plate the cells on appropriate medium and antibiotic. For cells transformed with the control pUC19 DNA, plate on LB agar containing 100 µg/mL of ampicillin. The remaining cell outgrowth can be stored at 4 °C in the event additional cell dilutions are plated.

6. Transformation of TransforMax EPI300 Chemically Competent *E. coli*

This procedure was written for the transformation of 50 µL of TransforMax EPI300 Chemically Competent *E. coli*. The procedure can be scaled down as needed.

1. Prepare 250 µL of SOC medium (**do not** include antibiotic in the medium) for each transformation to be performed. Maintain the medium at room temperature.
2. Chill 1.5 mL microcentrifuge tubes or microplate at 4 °C and heat a water bath or other temperature-controlled apparatus to 42 °C.
3. Thaw TransforMax EPI300 Chemically Competent *E. coli* cells on ice. Mix by gentle tapping. Use the cells immediately.
NOTE: *Refreezing chemically competent cells will result in a greatly reduced transformation efficiency.*
4. Transfer 1-5 µL of DNA and 50 µL of cells to a pre-chilled, microcentrifuge tube or into the wells of a microplate. Cover the cells and incubate on ice for 30 minutes.
5. Transfer the tubes or plate to 42 °C and heat shock for 30 seconds.
6. Transfer the cells back to ice and cool for 2 minutes.
7. Add 250 µL of the SOC medium to each tube.
8. Recover the cells by incubating at 37 °C for 60 minutes with shaking at 220-230 rpm.
9. Plate the cells on the appropriate media and antibiotic.

7. Induction of CopyControl fosmid clones

CopyControl Fosmid clones grown in TransforMax EPI300 cells can be amplified to 10-50 copies per cell. The induction process can be done in any culture volume desired depending on the needs of the user. Generally, a 1 mL induced culture will provide a sufficient amount of DNA for most applications

Manual

TransforMax EPI300 Electrocompetent and Chemically Competent *E. coli*

including sequencing and fingerprinting. Here we provide the procedure for amplifying the clones in 1 ml, 5 ml and 50 ml cultures.

Growth media for CopyControl fosmid clones

LB + chloramphenicol (12.5 µg/mL)

1. Add 5 mL of the appropriate Growth Media to 15 mL tubes for each fosmid or PCR clone that will be induced to high copy number.
2. Individually inoculate the media with a small portion of the desired fosmid or PCR clones grown on an overnight plate.
3. Grow the cultures overnight at 37 °C with shaking. These cultures will be used as inocula for the copy number amplification procedure.
4. From the table below, combine the appropriate volumes of fresh Growth media, the overnight culture and the CopyControl Induction Solution for the desired volume of induction culture. Aeration of the induction cultures is critical. Therefore, to maximise the surface area of the culture solution in the tube, perform the induction in the largest volume tubes that reasonably meets your needs and resources. For example, induce clones to high copy number in 1 mL of culture, using 1.5 mL tubes or larger, 5 mL cultures in 15 mL tubes and 50 mL cultures in 125 mL flasks.

Total volume of clone induction culture	Volume of fresh LB + chloramphenicol (12.5 µg/mL)	Volume of overnight 5 ml culture	Volume of 1000X CopyControl Induction Solution [§]
1 mL	800 µL	200 µL	1 µL
5 mL	4.5 mL	500 µL	5 µL
50 mL	45 mL	5 mL	50 µL

[§] Mix thoroughly after thawing.

5. Vigorously shake the tubes at 37 °C for 5 hours. Aeration is critical! Shake the tubes in a manner that will maximise aeration of the cultures (for example 1.5 mL tubes can be taped horizontally to the shaking table).
6. Centrifuge the cells and purify the DNA by your standard lab methods.

8. References

1. Hanahan D, 1983, *J. Mol. Biol.*, 166, 557.

9. Further support

If you require any further support, please do not hesitate to contact our Technical Support Team: techsupport@lgcgroup.com.

* TransforMax™ EPI300™ *E. coli* and related CopyControl™ Products are covered by U.S. Patent No. 5,874,259 licensed to LGC, Biosearch Technologies and/or by other patents pending and assigned to Biosearch Technologies for specific CopyControl vectors and/or cell lines. By purchasing CopyControl systems or vectors, the purchaser receives the right to use the product purchased from Biosearch Technologies or clones obtained therefrom solely for life science research. Purchase of these products does not grant rights to: (1) offer products, components of products, or any derivatives thereof for resale; or (2) to distribute or transfer the products, components of products, or derivatives thereof to third parties.



Integrated tools. Accelerated science.

   @LGCBiosearch | biosearchtech.com

All trademarks and registered trademarks mentioned herein are the property of their respective owners. All other trademarks and registered trademarks are the property of LGC and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details. No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopying, recording or any retrieval system, without the written permission of the copyright holder. © LGC Limited, 2021. All rights reserved. GEN/1018/EK/1221

For Research Use Only. Not for use in diagnostic procedures.



BIOSEARCH™
TECHNOLOGIES

GENOMIC ANALYSIS BY LGC