

# MC1061 F- Electrocompetent Cells



IMPORTANT!

**-80°C Storage Required**

Immediately Upon Receipt

Lucigen Corporation 2905 Parmenter St, Middleton, WI 53562 USA  
Toll Free: (888) 575-9695 | (608) 831-9011 | FAX: (608) 831-9012  
[lucigen@lucigen.com](mailto:lucigen@lucigen.com) [www.lucigen.com](http://www.lucigen.com)

FOR RESEARCH USE ONLY. NOT FOR HUMAN OR DIAGNOSTIC USE  
MA137 Rev. A, DCO No. 12-531

# MC1061 F- Electrocompetent Cells

## Contents

Components & Storage Conditions.....	2
MC1061 F- Electrocompetent Cells .....	3
MC1061 F- Electrocompetent Cells Transformation Protocol .....	3
Related Products.....	4

## Components & Storage Conditions

MC1061 F- Electrocompetent Cells are packed in Lucigen's DUO format for your convenience:

**DUO:** 50 µL per vial, enough for two 25-µL reactions

The cells are shipped on dry ice in one container, along with Recovery Medium and supercoiled control pUC19 DNA at 10 pg/µl. Please refer to Table 1 for a complete listing of MC1061 F- Electrocompetent Cells, efficiencies, and catalog numbers.

**Competent Cells require storage at  $-80^{\circ}\text{C}$ .**



**Table 1: MC1061 F- Competent Cells available from Lucigen**

STRAIN	Format	Efficiency (cfu/µg pUC19)	Transformations		Catalog #	Storage
MC1061 F- Electrocompetent Cells	DUO	$> 4 \times 10^{10}$	12 24	( 6 x 50 µL ) ( 12 x 50 µL )	60514-1 60514-2	<b><math>-80^{\circ}\text{C}</math></b>
Recovery Medium	-		12 24	( 1 x 12 mL ) ( 2 x 12 mL ) ( 8 x 12 mL )	---- ---- 80026-1	<b><math>-80</math> to <math>+20^{\circ}\text{C}</math></b>
*Supercoiled pUC19 DNA (10 pg/µL)	-			( 1 x 20 µL )	----	<b><math>-80</math> to <math>-20^{\circ}\text{C}</math></b>

\*Provided as a positive control for transformation.

# MC1061 F- Electrocompetent Cells

## MC1061 F- Competent Cells

MC1061 F- is a non amber-suppressing strain that is commonly used for general cloning and Phage display because of its ability to reach very high transformation efficiencies. It is identical to Lucigen strain SS320, except that it lacks the F' episome that is required for infection by filamentous phage such as M13. The strain can be used for plasmid propagation and phage display, but cannot be used for re-infection.

### Transformation Efficiency:

Electrocompetent Cells:  $\geq 4 \times 10^{10}$  cfu/ $\mu$ g

### Genotype

*araD139 Δ (araA-leu)7697 Δ(lac)X74 galK16 galE15(GalS) lambda- e14- mcrA0 relA1 rpsL150 (Str<sup>R</sup>) spoT1 mcrB1 hsdR2*

## Usage Guidelines

### Preparation for Transformation

Transformation is carried out in a 0.1 cm gap cuvette using 25  $\mu$ L of MC1061 F- Electrocompetent Cells. Optimal settings for electroporation are listed in the table below. Typical time constants are 3.5 to 4.5 msec.

Optimal Setting	Alternate Settings (~ 20-50% lower efficiencies)	For 2.0 mm cuvettes
1.0 mm cuvette 10 $\mu$ F 600 Ohms 1800 Volts	1.0 mm cuvette 25 $\mu$ F 200 Ohms 1400 – 1600 Volts	2.0 mm cuvette 25 $\mu$ F 750 Ohms 2400 Volts

### Suggested Electroporation Systems:

Bio-Rad Micro Pulser #165-2100; Bio-Rad E. coli Pulser #165-2102; Bio-Rad Gene Pulser II #165-2105; BTX ECM630 Electroporation System; Eppendorf Electroporator 2510.

Optional transformation control reactions include electroporation with 1  $\mu$ L (10 pg) of supercoiled pUC19 DNA.

To ensure successful transformation results, the following precautions must be taken:

- Ligation reactions performed with Lucigen's CloneDirect™ Ligation Buffer (included with Lucigen's Cloning or Ligation Kits) must be heat killed at 70°C for 15 minutes before transformation. The CloneDirectligation reaction can be used directly after heat inactivation, without purification of the ligation products.
- DNA samples in other buffers must be purified and dissolved in water or a buffer with low ionic strength, such as TE. The presence of salt in the electroporation sample leads to arcing at high voltage, resulting in loss of the cells and DNA.
- Microcentrifuge tubes and electroporation cuvettes must be thoroughly pre-chilled on ice before use. Successful results can be obtained with cuvettes from many sources including BTX (Model 610), Eppendorf (Cat. #940001005), and BioRad (Cat. #165-2089).
- The cells must be completely thawed **on ice** before use.

# MC1061 F- Electrocompetent Cells

---

- For highest transformation efficiency, use the provided Recovery Medium to resuspend the cells after electroporation. Use of SOC or other media will result in lower transformation efficiencies.
- Prepare nutrient agar plus appropriate antibiotic.

## Transformation Protocol

1. Have Recovery Medium and 17 mm x 100 mm sterile culture tubes readily available at room temperature (one tube for each transformation reaction). Transformation efficiency may decrease with the use of SOC or other media.
2. Place electroporation cuvettes (0.1 cm gap) and microcentrifuge tubes on ice (one cuvette and one microfuge tube for each transformation reaction).
3. Remove cells from the -80°C freezer and place on wet ice until they thaw **completely** (10-20 minutes).
4. When cells are thawed, mix them by tapping gently. Aliquot 25 µL of cells to the chilled microcentrifuge tubes on ice.
5. If using ligation buffer from any Lucigen Cloning Kit, add 1 µL of the heat-denatured ligation reaction to the 25 µL of cells on ice. Failure to heat-inactivate the ligation reaction will prevent transformation. Stir briefly with pipet tip; **do not** pipet up and down to mix, which can introduce air bubbles and warm the cells. Use of more than 2 µL of ligation mix may cause electrical arcing during electroporation.

For ligation reactions using other commercial kits, please refer to the manufacturer's instructions.

6. Carefully pipet 25 µL of the cell/DNA mixture into a chilled electroporation cuvette without introducing bubbles. Quickly flick the cuvette downward with your wrist to deposit the cells across the bottom of the well. Electroporate according to the conditions recommended above.
7. Within 10 seconds of the pulse, add 975 µL of Recovery Medium to the cuvette and pipet up and down three times to resuspend the cells. Transfer the cells and Recovery Medium to a culture tube.
8. Place the tube in a shaking incubator at 250 rpm for 1 hour at 37°C.
9. Spread up to 100 µl of transformed cells on nutrient agar plates containing the appropriate antibiotic.
10. Incubate the plates overnight at 37°C.
11. Transformed clones can be further grown in any rich culture medium.

## Related Lucigen Products

- NxGen™ Terminal Transferase
- NxGen™ T7 RNA Polymerase
- DNATerminator® End Repair Kit
- PCRTerminator® End Repair Kit
- UltraClone™ DNA Ligation & Transformation Kits
- CloneDirect™ Rapid Ligation Kit
- PCR-SMART™ Cloning Kits
- *E. coli*® EXPRESS Electrocompetent Cells
- OverExpress™ Competent Cells

**Notice of Limited Label License, Copyright, Patents, Warranties, Disclaimers, and Trademarks**

# MC1061 F- Electrocompetent Cells

---

Copyright© 2012 by Lucigen Corp. All rights reserved. Lucigen, BigEasy, ClonePlex, CloneSmart, CopyRight, DNATerminator, E. cloni, EconoTaq, Expresso, HI-Control, pETite, PCR Terminator, pJAZZ, pSMART, PyroPhage, ChimeraFree, CloneDirect, cSMART, Expresso, eLucidations, GapFree, NxGen, NxSeq, PCR SMART, pEZ, pEZSeq, pGC, pRham, pLEXX-AK, pRANGER, Replicator, Taq98, TaqSelect, The Cloning Company, The Molecular Cloning Company, Transformance, TSA, PyroScript, Endura and ClonID are registered trademarks of Lucigen Corporation.

Lucigen's products are sold for research use only and are not to be used in humans or for medical diagnostics. Lucigen's liability with respect to any of its products is limited to the replacement of the product. No other warranties of any kind, expressed or implied, including without limitation, any implied fitness for any particular use, are provided by Lucigen. Lucigen is not liable for any direct, indirect, incidental or consequential damages arising out of or in connection with the use or inability to use any of its products.

If the purchaser is not willing to accept these use limitations, Lucigen Corporation is willing to accept return of the product for a full refund. For information on obtaining a license, contact Lucigen Corporation, 2905 Parmenter St., Middleton, WI 53562. Email: [Lucigen@lucigen.com](mailto:Lucigen@lucigen.com). Phone: 608-831-9011. Fax 608-831-9012.

## **Warranty**

Lucigen warrants that its products will conform to the standards stated in its product specification sheets in effect at the time of shipment. Lucigen will replace, free of charge, any product that does not conform to the specifications. This warranty limits Lucigen's liability only to the replacement of the product. THIS WARRANTY IS EXCLUSIVE, AND LUCIGEN MAKES NO OTHER WARRANTY, EXPRESS OR IMPLIED, INCLUDING WITHOUT LIMITATION, ANY IMPLIED WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE. The stated express warranties, and the remedy provided for breach thereof, are in lieu of all other liability or obligations of Lucigen for any damages whatsoever arising out of or in connection with the delivery, use, performance, or the inability to use any of its products. IN NO EVENT SHALL LUCIGEN BE LIABLE UNDER ANY LEGAL THEORY (INCLUDING BUT NOT LIMITED TO CONTRACT, NEGLIGENCE, STRICT LIABILITY IN TORT, OR WARRANTY OF ANY KIND) FOR ANY INDIRECT, SPECIAL, INCIDENTAL, CONSEQUENTIAL, OR EXEMPLARY DAMAGES (INCLUDING BUT NOT LIMITED TO LOST PROFITS). Without limiting the effect of the preceding sentence, Lucigen's maximum liability, if any, shall not exceed the purchase price paid by buyer for the product.