

Manual

Bst DNA Polymerase, Exonuclease Minus 50,000 units/mL

For Research Use Only. Not for use in diagnostic procedures.

IMPORTANT
-20 °C storage required
immediately upon receipt

Manual

Bst DNA Polymerase, Exonuclease Minus 50,000 units/mL

1. Product description

Bst DNA Polymerase, Exonuclease Minus, 50,000 units/mL.

2. Product specifications

Storage buffer: 10 mM Tris-HCl, pH 7.5, 50 mM KCl, 1 mM DTT, 0.1 mM EDTA, 0.1% Triton® X-100 (Rohm & Haas) and 50% Glycerol.

Stability: Bst DNA Polymerase, Exonuclease Minus is stable for one year from the date received if stored at -20 °C.

Recommended reaction conditions: 50 units Bst DNA Polymerase, Exonuclease Minus; 1X DNA Polymerase Buffer B containing 20 mM Tris-HCl pH 8.8, 10 mM (NH₄)₂SO₄, 10 mM KCl, 2 mM MgSO₄ and 0.1 % Triton X-100.

Activity determination: One unit catalyses the incorporation of 10 nmol of dNTP into acid-insoluble material in 30 minutes at 65 °C in 20 mM Tris-HCl pH 8.8, 10 mM (NH₄)₂SO₄, 10 mM KCl, 2 mM MgSO₄, 0.1 % Triton X-100, 30 nM M13mp18 ssDNA, 70 nM M13 sequencing primer(-47) 24 mer, 200 µM dGTP, dATP, dTTP, dCTP (a mix of unlabeled and [³³P]dCTP) and 0.1 mg/ml BSA.

Absence of endonuclease or nicking activity: Incubation of 50 units of Bst DNA Polymerase, Exonuclease Minus with 1 µg of pUC19 DNA for 16 - 18 hours at 37 °C resulted in no smearing of bands detected by agarose gel electrophoresis.

Absence of exonuclease activity: Incubation of 50 units of Bst DNA Polymerase, Exonuclease Minus with 1 µg of *Hind*III-cut lambda DNA for 16 hours at 37 °C and 65 °C resulted in no smearing of bands on agarose gels. Single stranded and double stranded exonuclease activities were tested by incubating 10 µL of enzyme at 50 units/µL with radiolabeled DNA substrate for one hour at 37 °C and 65 °C, resulting in less than 0.1% release of TCA-soluble counts.

Purity: >90% pure by SDS PAGE. No detectable DNA contamination. 10 µL of enzyme at 50 units/µL of the sample was tested for *E. coli* genomic DNA contamination by PCR amplifying with the *E. coli* 16S ribosomal primers.

Heat inactivation: 80 °C for 20 minutes.

3. Product designations and kit components


Product	Concentration	Kit size	Catalogue number	Reagent description	Part number	Volume
Bst DNA Polymerase, Exonuclease Minus 50,000 units/ml	50,000 units/mL	10,000 units	30028-1	Bst DNA Polymerase, Exonuclease Minus (50 units/µL)	F93636-1	200 µL
				10X DNA Polymerase Buffer B	F98637-1	4.8 mL

Manual

Bst DNA Polymerase, Exonuclease Minus 50,000 units/mL

4. Storage conditions

Store all kits and components at -20 °C

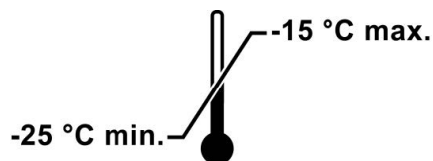

5. Applications

1. DNA sequencing through high GC regions^{1,2}
2. Rapid Sequencing from nanogram amounts of DNA template³

6. References

1. Griffin H and Griffin A 1994 *PCR Technology*, 228-229.
2. McClary J *et al.* 1991 *J. DNA Sequencing and Mapping*, 1, 173-180
3. Mead DA *et al.* 1991 *Biotechniques*, 11, 76-87.

7. Technical support

If you require any further support, please do not hesitate to contact our Technical Support Team:

techsupport@lgcgroup.com.

Legal and warranty

LGC Biosearch Technologies™ does not encourage or support the unauthorized or unlicensed use of patented nucleic acid amplification processes. It is the sole responsibility of the buyer to ensure that use of the product does not infringe the patent rights of third parties. If the purchaser is not willing to accept these use limitations, Biosearch Technologies is willing to accept return of the product for a full refund.

Biosearch Technologies warrants that its products will conform to the standards stated in its product specification sheets in effect at the time of shipment. Biosearch Technologies will replace, free of charge, any product that does not conform to the specifications. This warranty limits Biosearch Technologies' liability only to the replacement of the product. THIS WARRANTY IS EXCLUSIVE, AND BIOSEARCH TECHNOLOGIES MAKES NO OTHER WARRANTY, EXPRESS OR IMPLIED, INCLUDING WITHOUT LIMITATION, ANY IMPLIED WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE. The stated express warranties, and the remedy provided for breach thereof, are in lieu of all other liability or obligations of Biosearch Technologies for any damages whatsoever arising out of or in connection with the delivery, use, performance, or the inability to use any of its products. IN NO EVENT SHALL BIOSEARCH TECHNOLOGIES BE LIABLE UNDER ANY LEGAL THEORY (INCLUDING BUT NOT LIMITED TO CONTRACT, NEGLIGENCE, STRICT LIABILITY IN TORT, OR WARRANTY OF ANY KIND) FOR ANY INDIRECT, SPECIAL, INCIDENTAL, CONSEQUENTIAL, OR EXEMPLARY DAMAGES (INCLUDING BUT NOT LIMITED TO LOST PROFITS).



   @LGCBiosearch | biosearchtech.com

All trademarks and registered trademarks mentioned herein are the property of their respective owners. All other trademarks and registered trademarks are the property of LGC and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details. No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopying, recording or any retrieval system, without the written permission of the copyright holder. © LGC Limited, 2023. All rights reserved. GEN/954/EK/0123

For Research Use Only. Not for use in diagnostic procedures.



BIOSEARCH™
TECHNOLOGIES

GENOMIC ANALYSIS BY LGC