

Product insert

Protease K solution 20 mg/mL

For Research Use Only. Not for use in diagnostic procedures.

Product overview

Protease K solution is provided in solution at 20 mg/mL. Enzymatic activity is ≥ 600 U/mL. The solution is supplied at pH 8.0.

Protease K is stable over a wide pH range (4.0-12.5), with optimal pH range of 7.5-8.0.

Protease K is most active at 55-60 °C, and functions optimally at 56 °C. A decrease in activity may be observed at temperatures above 60 °C. A suggested incubation would be 30 minutes at 56-60 °C. When working in plates, we recommend incubating at 60 °C as heat transfer in plates is not optimal and it can take time until the sample(s) themselves reach 60 °C.

Please note that RNase and DNase can be digested by Protease K when added simultaneously. We recommend adding RNase to a sample first, incubating as required, before the Protease K is added and a subsequent incubation performed. Exact incubation times will need to be validated for your specific application. The enzyme can be completely inactivated through incubation at 95 °C for 10 minutes.

Storage

Protease K solution should be stored at room temperature and is stable for 1 year. Please refer to product label for exact expiry date. The product can be shipped at room temperature. It is a very heat-stable and robust enzyme, and long-term storage of our protease solution at higher than ambient temperatures has shown no reduction in activity.

Ordering information

Cat no.	Description	Volume
NAP30-002-01	Protease K solution 20 mg/mL	1 mL
NAP30-002-02	Protease K solution 20 mg/mL	10 mL
NAP30-002-03	Protease K solution 20 mg/mL	100 mL
NAP30-002-04	Protease K solution 20 mg/mL	250 mL
NAP30-002-05	Protease K solution 20 mg/mL	500 mL
NAP30-002-06	Protease K solution 20 mg/mL	1000 mL

Safety information

The SDS for Protease K solution can be accessed via our web shop: shop.biosearchtech.com.




 @LGCBiosearch | biosearchtech.com

All trademarks and registered trademarks mentioned herein are the property of their respective owners. All other trademarks and registered trademarks are the property of LGC and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details. No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopying, recording or any retrieval system, without the written permission of the copyright holder. © LGC Limited, 2023. All rights reserved. GEN/1074/FS/0123